

ADDITIONAL PERISTALTIC PUMP

CAB-2

DOUBLING THE FLOW DURING
DISPENSING

ACCESSORY



Additional peristaltic pump

APPLICATION

- + This option is perfect for applications that require the dispensing of small to medium volumes, such as the production of test tubes in microbiology, the filling of Petri dishes, and other similar processes.



INCLUDED ITEMS

- 1 silicone tube, 2m long and $\varnothing 6,4$ (inside) x $\varnothing 9,6$ mm (outside).
- 1 complementary head.
- 1 Y-shaped tube.

BENEFITS

- Increase in productivity.
- Easy installation at any time.

OPERATION

To activate each dispensation, there are two available mechanisms:

- Mode 1: Manual dispensing using a foot pedal.
- Mode 2: Semi-automatic dispensing with delay.

Depending on the combination of the chosen tube size and the number of installed peristaltic pumps, the dispensing speed in this mode ranges between 7 and 33mL/s:

Other tube diameters available.



DISPENSING LINES

Ref. TUB-DOSIF

Available diameters: 3,2; 4; 4,8; 6,4 and 8mm.

Dispensing performance	Ø tube (mm)	Dispensing speed
Dispensing using a single peristaltic pump	3,2	7mL/s
	4	9mL/s
	4,8	11mL/s
	6,4	15mL/s
	8	20mL/s
Dispensing using two peristaltic pumps	3,2	12mL/s
	4	15mL/s
	4,8	18mL/s
	6,4	25mL/s
	8	33mL/s




Additional peristaltic pump


INSTALLATION

It can be installed at any time.


1. Remove the head of the equipment.

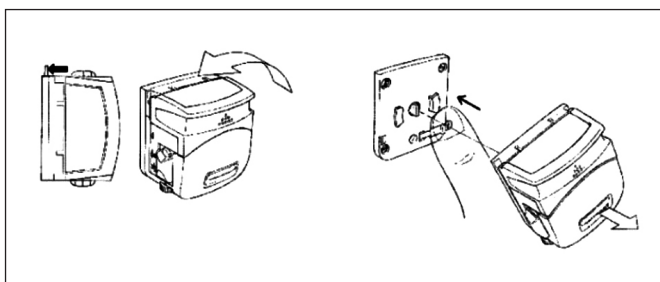

2. Place the complementary head CAB-2 in its place.

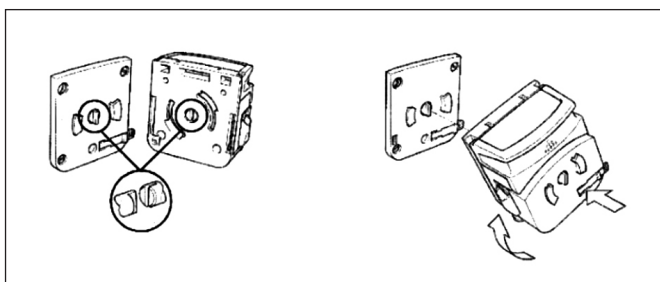

3. Fit the equipment's head into the complementary head.

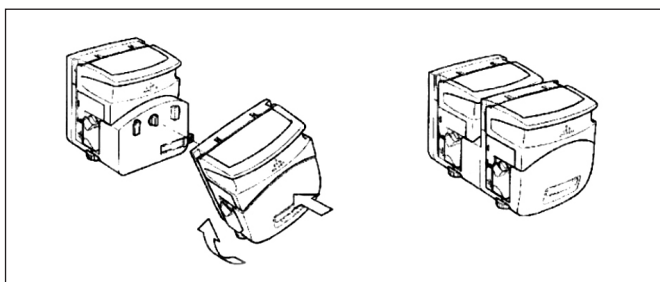

REV 03.2025

